

NEULOG BAROMETER LOGGER SENSOR GUIDE



NeuLog barometer logger sensor NUL-221

The NeuLog barometer sensor can be used for any science experiment or activity which requires an accurate atmospheric pressure measurement such as in the fields of Physics, Biology, Physiology, Exercise Science, Environmental Science, Earth Science, Weather Science, Botany, etc.

The barometer sensor comes pre-calibrated in the factory, and never requires recalibration.

Just a few of the thousands of possible experiments that can be done with the NUL-221 sensors are: demonstrating how pressure is related to altitude, exercise studies at different altitudes, effects of pressure change on plant growth, weather experiments, and many more.

The pressure sensor can collect data using four different units of measure:

- Kilopascal (kPa): The SI unit of pressure.
- Atmospheres (atm): A pressure unit equal to one earth atmosphere at sea level (101.325 kPa)
- Inches of mercury (in.Hg): A non-SI unit of atmospheric pressure
- Millimeters of mercury (mmHg): A non-SI unit of atmospheric pressure

Quick start procedure:

PC or Mac Computer

Materials needed:

- NUL-221 Barometer Sensor
- USB-200 USB Module
- A USB to mini USB cable (which comes with the USB-200)

Your barometer sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via a USB to mini-USB cable. Please note that you cannot plug the barometer sensor directly into the computer.

Resident PC software and browser based software can be downloaded for free at www.NeuLog.com/download as well as a full software user guide. **Note:** Make sure not to download and install both types of software, they will conflict on the computer.

Procedure:

1. Install the NeuLog software
2. Connect the USB-200 module to the PC or Mac
3. Connect the barometer sensor to the USB-200 module (they directly plug together). Please note there is no calibration required for this sensor.
4. Open the NeuLog software.
5. Once a barometer sensor logo appears on the left side of the screen the probe has been automatically identified and you can begin experimentation.
6. If the barometer sensor is not automatically identified then click the "Search for sensors" icon to find the sensor.
7. Select the "On-line experiment" button; this will open a graph below.
8. Click on the "Module setup" button located on the barometer sensor icon in the module window to change the sensor settings if need be.
9. Click on the experiment set up button to change the experiment settings if need be (experiment duration for example).
10. The barometer sensor will give a live reading in the box to the left of the screen while plugged in.
11. To run an experiment and collect data click "Run experiment".
12. To end data collection early, click "Stop experiment".

Tablet, smart phone device

Materials needed:

- NUL-221 Barometer Sensor
- WIFI-201 WIFI module
- BAT-200 Battery

Your barometer sensor needs to be connected to a WIFI-201 module. The WIFI-201 module will create a closed NeuLog wifi network which will stream the NeuLog data to a device of your choosing. Once your device is wirelessly connected to the NeuLog network you can run experiments and collect data through a browser of your choosing.

Procedure:

1. Connect the barometer sensor directly to the left side of a WIFI-201 module (no wires required).
2. Connect a BAT-200 module to the right side of the WIFI-201 module.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to micro USB charger (such as a typical cell phone charger).

NEULOG BAROMETER LOGGER SENSOR GUIDE



The WIFI-201 module will run for 15-75 minutes (depending on the sensor) without being plugged in.

- For further WIFI-201 instructions or the WIFI-201 quick start guide please visit www.NeuLog.com/download.
- The WIFI-201 indicator lights will flash; take no action until the LED to the far left turns blue.
- Take your tablet or smart phone and go to the Wi-Fi settings and select the NeuLog network which matches the WIFI module ID found on the back of the WIFI-201 device. (NeuLogXXXX for example)
- Give your device 1-2 minutes to connect to the WIFI-201 network.
- Once the device is connected go to your browser and type in the website wif201.com into the URL bar, then wait for 30-60 seconds.
- You will see a "Control mode" icon in the browser, click on this icon.
- The browser will then load a new screen and begin to auto detect the sensors, this can take a few minutes.
- If the browser does not auto detect, select "Search for sensors"
- Once the sensor is found you will see an icon on the left side of the screen for the barometer sensor. The icon will display data in real time.
- To set the barometer sensor's settings click on "Module setup" button located on the barometer sensor icon on the left side of your screen.
- To change the experimental settings click the "Experiment setup" button (experiment duration for example).
- To run an experiment and collect data click "Run experiment".

Operation with NeuLog viewer

Materials needed:

- NUL-221 Barometer Sensor
- VIEW-101 Viewing Module
- BAT-200 Battery

Procedure:

- Connect the barometer sensor to the left side of the viewer.
- Connect the BAT-200 module to the right side of the viewer.
- The VIEW-101 will auto detect the sensor and you will see it appear on the left side of the screen.
- Once the sensor appears it will be monitoring data in real time.

- To run an experiment and collect data click the run button (little green person).

Off-line experiments

(Off-line experiments are for when you do not have a sensor connected directly to a computer, tablet, smartphone, or NeuLog viewer).

Materials needed:

- NUL-221 Barometer Sensor
- BAT-200 Battery

Materials needed to configure your offline experiment:

- USB-200 USB Module or WIFI-201 WIFI module or VIEW-101 Viewing Module
- A USB to mini USB cable (which comes with the USB-200)

Procedure:

- Connect the barometer sensor directly to the left side of a charged BAT-200 module.
- When ready to collect data press the "Start/Stop" button on the barometer sensor, a red light will turn on for the duration of the experiment.
- To change the experiment settings first connect to your chosen device (PC, Mac, tablet, smart device) and select the "Module setup" button. (For more information on how to change the experiment settings view quick start procedure section for your chosen device in this document.)
- After the experiment has concluded (when the "Start/Stop" button is pressed again or the data collection period ends) connect the sensor to your chosen device. (full instructions in each device's quick start procedure section)
- Open NeuLog software.
- Click the "Off-line experiment" button.
- Click the "Load data from sensors" button.
- Select which experimental data to upload. (5 experimental runs can be stored at one time on the NUL-221 barometer sensor)

NEULOG BAROMETER LOGGER SENSOR GUIDE



Restoring sensor's factory default settings:

Procedure:

1. Connect the USB-200 to a PC, Mac, tablet, or smart device.
2. Connect the barometer sensor to a USB-200 module (they directly plug together). Please note there is no calibration required for this sensor.
3. Open the NeuLog software.
4. Click the "Tools" icon.
5. Click "Restore sensor's factory defaults" icon.
6. If prompted to clear the graph, click either the "Off-line experiment" or "On-line experiment" (whichever you are using).
7. Click "Clear experiment results".
8. Resume from step 4.

Included with sensor:

- NeuLog NUL-221 Barometer Sensor Guide (this document).

Sensor specifications:

Range and operation modes	ADC resolution	Resolution	Max sample rate (S/sec)
80 to 106 kPa	15 bit	0.1 kPa	100
0.8 to 1.00 atm		0.01 atm	
23.60 to 31.30 in Hg		0.01 in Hg	
600 to 795 mm Hg		0.7 mm Hg	
-380 to 1950 m		0.1 m	

Sensor features:

- Fully digital data.
- Rugged plastic ergonomic case.
- Push button switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).
- Pre-calibrated sensing equipment.
- **Note:** NeuLog products are intended for educational use.

Videos and experiment examples:

- Videos, literature and other probes can be found at www.NeuLog.com
- <http://www.NeuLog.com/Products/barometerloggerSensor.aspx>
- **Air Pressure and Altitude -**
<http://www.neuLog.com/Portals/3/Users/Experiment E-2 Air Pressure and Altitude Ver 3.0.2.pdf>

Technical background:

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (°C, °F, Lux, %, ppm, for example).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

Inside the barometer sensor's plastic body there is a flexible membrane which bends with atmospheric pressure differences. On one side of the membrane there is a vacuum which is compared against the other side of the membrane, the atmosphere. An electrical charge is generated based on the amount of flexing in the membrane. Temperature is taken into account using internal circuitry during the pressure calculation as to normalize the data for any thermal fluctuations.

Maintenance and storage:

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the barometer sensor's body.
- After use, gently wipe away any foreign material from the barometer sensor.
- Store in a box at room temperature out of direct sunlight.

Warranty:

We promise to deliver our sensor free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your sensor will be repaired or replaced.

NEULOG BAROMETER LOGGER SENSOR GUIDE



Thank you for using NeuLog!



Flexible, simple, fast, forward thinking.

W: www.neulog.com

E: info@neulog.com

A: 850 St Paul Street, Suite 15, Rochester, NY 14605

P: 1.866.553.8536

V04302013