

NEULOG ELECTROCARDIOGRAM LOGGER SENSOR GUIDE



NeuLog electrocardiogram Logger Sensor NUL-218

The NeuLog electrocardiogram sensor can be used for any science experiment or activity that can utilize heartbeat data such as in the fields of: Biology, Physiology, Human medicine, Psychology, etc.

An electrocardiogram is a very commonly used medical device which measures and graphs the heart's electrical activity. Every single heartbeat is a very complex event which involves many anatomical locations undergoing changes in their electrical charges. A depolarizing cascade event originates in the Sinoatrial Node, which is often referred to as the natural pacemaker, and travels throughout the rest of the heart.

The sensor comes pre-calibrated so you can start experimentation right out of the box using any of the following guides.

Just a few of the thousands of possible experimental subjects that can be done with the NUL-218 sensors are: electrical impulses, heart rate monitoring, human physiology, muscle physiology, heart rate changes to stimuli, and many more.

The ECG sensor uses the following unit of measure:

- Arbitrary analog units (Arb): An arbitrary unit to demonstrate waves, frequencies, and periods. Therefore what you will see on the Y-axis is signal intensity, while the X-axis shows time.


Electrocardiograms:

The heart undergoes a very complex process of electrical activity; this is the basis for how electrocardiograms work. As the heart's muscles contract, the cells are electrically depolarized which causes a cascade of depolarization to flow across the heart – contracting the heart's muscle fibers.

During a heartbeat a very complex series of events happens which begins with the cells resting potential at roughly -90mV (millivolts). The heart's Sinoatrial Node begins the cascade by depolarizing – creating an electrical current. The electrical signal

travels to the right atrium and almost immediately to the left atrium. Because the atria and the ventricles are insulated from each other by a special type of cell which acts as an electrical insulator, the signal is picked up in another cluster of cells called the Atrioventricular Node located in the right atria. The signal is carried by a special type of fibers - the Bundle of His into the ventricles. Purkinje Fibers rapidly spread the electrical depolarization cascade throughout the ventricles. The entire event takes some time to complete. The graph is a visual demonstration of the heart's relative charge at any given point during heartbeats.

The electrical current is measured on the skin using conducting leads and converted into a graphic that is very standard and reproducible.

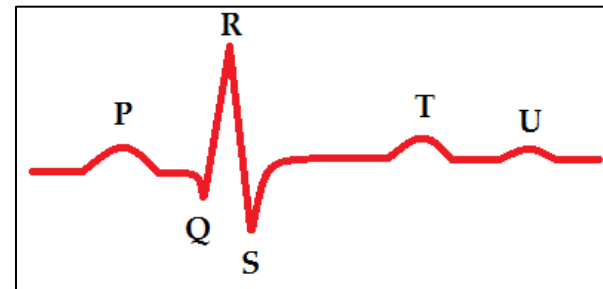

- P:** Contraction of atria
- Q:** Initial depolarization of the ventricle
- R:** Early depolarization of the ventricles
- S:** Late depolarization of the ventricles
- T:** Repolarization of the ventricles
- U:** Often hidden by the T wave – this wave has an unknown origin

This graph represents one complete heart beat from depolarization of the Sinoatrial Node to repolarization of the ventricles.

Connecting to the ECG sensor:

The NeuLog electrocardiogram sensor uses three reusable clip leads for easy connection to the arms. Each pair of leads (red + yellow, red + black, black + yellow) measure the heart's electrical fluctuations and are all compared against one another.

NEULOG ELECTROCARDIOGRAM LOGGER SENSOR GUIDE



Leads:

- Black: Ground
- Red: Positive
- Yellow: Negative

How to connect:

1. Connect the NeuLog ECG sensor to a smart device using one of the following guides. (PC or Mac/Tablet/Viewer)
2. Place the red clip on the subject's left wrist with the metal connector on the underside (palm side) of the arm.
3. Place the yellow clip on the subject's right wrist with the metal connector on the underside (palm side) of the arm.
4. Place the black clip on the subject's left forearm or elbow.
5. Connect the three alligator clips (red, yellow, black) directly to their corresponding arm clip's (red, yellow, black) metal piece that forms a loop.
6. You are now connected.

Note: There are other configurations possible, this is just one possibility. If the R wave shows a downward slope instead of upward, your connection's polarization is reversed though the absolute results are still the same.

Quick start procedure:

PC or Mac Computer

Materials needed:

- NUL-218 ECG Sensor
- 3 Neulog Reusable Clamp Leads
- USB-200 USB Module
- A USB to mini USB cable (which comes with the USB-200)

Your ECG sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via a USB to mini-USB cable. Please note that you cannot plug the ECG sensor directly into the computer.

Resident PC software and browser based software can be downloaded for free at www.NeuLog.com/download as well as a full software user guide. **Note:** Make sure not to download and install both types of software, they will conflict on the computer.

Procedure:

1. Install the NeuLog software
2. Connect the USB-200 module to the PC or Mac
3. Connect the ECG sensor to the USB-200 module (they directly plug together). Please note there is no calibration required for this sensor.
4. Open the NeuLog software.
5. Once an ECG sensor logo appears on the left side of the screen the probe has been automatically identified and you can begin experimentation.
6. If the ECG sensor is not automatically identified then click the "Search for sensors" icon to find the sensor.
7. Select the "On-line experiment" button; this will open a graph below.
8. Click on the "Module setup" button located on the ECG sensor icon in the module window to change the sensor settings if need be.
9. Click on the experiment set up button to change the experiment settings if need be (experiment duration for example).
10. The ECG sensor will give a live reading in the box to the left of the screen while plugged in.
11. To run an experiment and collect data click "Run experiment".
12. To end data collection early, click "Stop experiment".

Tablet, smart phone device

Materials needed:

- NUL-218 ECG Sensor
- 3 Neulog Reusable Clamp Leads
- WIFI-201 WIFI module
- BAT-200 Battery

Your ECG sensor needs to be connected to a WIFI-201 module. The WIFI-201 module will create a closed NeuLog wifi network which will stream the NeuLog data to a device of your choosing. Once your device is wirelessly connected to the NeuLog network you can run experiments and collect data through a browser of your choosing.

NEULOG ELECTROCARDIOGRAM LOGGER SENSOR GUIDE



Procedure:

1. Connect the ECG sensor directly to the left side of a WIFI-201 module (no wires required).
2. Connect a BAT-200 module to the right side of the WIFI-201 module.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to micro USB charger (such as a typical cell phone charger). The WIFI-201 module will run for 15-75 minutes (depending on the sensor) without being plugged in.
4. For further WIFI-201 instructions or the WIFI-201 quick start guide please visit www.NeuLog.com/download.
5. The WIFI-201 indicator lights will flash; take no action until the LED to the far left turns blue.
6. Take your tablet or smart phone and go to the Wi-Fi settings and select the NeuLog network which matches the WIFI module ID found on the back of the WIFI-201 device. (NeuLogXXXX for example)
7. Give your device 1-2 minutes to connect to the WIFI-201 network.
8. Once the device is connected go to your browser and type in the website wifi201.com into the URL bar, then wait for 30-60 seconds.
9. You will see a "Control mode" icon in the browser, click on this icon.
10. The browser will then load a new screen and begin to auto detect the sensors, this can take a few minutes.
11. If the browser does not auto detect, select "Search for sensors"
12. Once the sensor is found you will see an icon on the left side of the screen for the ECG sensor. The icon will display data in real time.
13. To set the ECG sensor's settings click on "Module setup" button located on the ECG sensor icon on the left side of your screen.
14. To change the experimental settings click the "Experiment setup" button (experiment duration for example).
15. To run an experiment and collect data click "Run experiment".

Operation with NeuLog viewer

Materials needed:

- NUL-218 ECG Sensor
- 3 Neulog Reusable Clamp Leads
- VIEW-101 Viewing Module

- BAT-200 Battery

Procedure:

1. Connect the ECG sensor to the left side of the viewer.
2. Connect the BAT-200 module to the right side of the viewer.
3. The VIEW-101 will auto detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data click the run button (little green person).

Off-line experiments

(Off-line experiments are for when you do not have a sensor connected directly to a computer, tablet, smartphone, or NeuLog viewer).\

Materials needed:

- NUL-218 ECG Sensor
- 3 Neulog Reusable Clamp Leads
- BAT-200 Battery

Materials needed to configure your offline experiment:

- USB-200 USB Module or WIFI-201 WIFI module or VIEW-101 Viewing Module
- A USB to mini USB cable (which comes with the USB-200)

Procedure:

1. Connect the ECG sensor directly to the left side of a charged BAT-200 module.
2. When ready to collect data press the "Start/Stop" button on the ECG sensor, a red light will turn on for the duration of the experiment.
3. To change the experiment settings first connect to your chosen device (PC, Mac, tablet, smart device) and select the "Module setup" button. (For more information on how to change the experiment settings view quick start procedure section for your chosen device in this document.)
4. After the experiment has concluded (when the "Start/Stop" button is pressed again or the data collection period ends) connect the sensor to your chosen device. (full instructions in each device's quick start procedure section)

NEULOG ELECTROCARDIOGRAM LOGGER SENSOR GUIDE



5. Open NeuLog software.
6. Click the "Off-line experiment" button.
7. Click the "Load data from sensors" button.
8. Select which experimental data to upload. (5 experimental runs can be stored at one time on the NUL-218 ECG sensor)

Restoring sensor's factory default settings:

Procedure:

1. Connect the USB-200 to a PC, Mac, tablet, or smart device.
2. Connect the ECG sensor to a USB-200 module (they directly plug together). Please note there is no calibration required for this sensor.
3. Open the NeuLog software.
4. Click the "Tools" icon.
5. Click "Restore sensor's factory defaults" icon.
6. If prompted to clear the graph, click either the "Off-line either" or "On-line experiment" (whichever you are using).
7. Click "Clear experiment results".
8. Resume from step 4.

Included with sensor:

- NeuLog NUL-218 ECG Sensor Guide (this document).
- Neulog ECG probe with three alligator clips for easy connection to clamp leads.
- 3 ECG clamp leads (red, yellow, black).

Sensor specifications

Sensor specifications	
Range and operation modes	0 to 4092 Arbitrary units
ADC resolution	12 bit
Resolution	1
Max sample rate (S/sec)	100

Sensor features:

- Fully digital data.
- Rugged plastic ergonomic case.
- Push button switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).
- ECG probe attached to sensor body by a durable rubber-coated wire with three alligator clip ends.
- Pre-calibrated sensing equipment.
- **Note:** NeuLog products are intended for educational use.

Videos and experiment examples:

- Videos, literature and other probes can be found at www.NeuLog.com
- <http://www.NeuLog.com/Products/ECGloggerSensor.aspx>
- **Electrocardiogram Measurement –**
<http://www.neulog.com/Portals/3/Users/Experiment B-40 Electrocardiogram Measurement Ver 3.0.2.pdf>

Technical background:

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (°C, °F, Lux, %, ppm, for example).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

ECG sensors detect minor changes in electrical voltage which originate in the heart and spread out across the body over a very short amount of time. Cells have a resting potential of about 0.9 mV. A heart beat originates from an electrical impulse in the Sinoatrial Node, often called the natural pacemaker. The electrical potential cascade spreads across the heart. Using three leads, a positive, a negative, and a ground the heart's electrical charge (which originated in the heart) is measured.

NEULOG ELECTROCARDIOGRAM LOGGER SENSOR GUIDE



Maintenance and storage:

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the ECG sensor's body.
- After use, gently wipe away any foreign material from the ECG sensor.
- Store in a box at room temperature out of direct sunlight.

Warranty:

We promise to deliver our sensor free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your sensor will be repaired or replaced.

Thank you for using NeuLog!



Flexible, simple, fast, forward thinking.

W: www.neulog.com

E: info@neulog.com

A: 850 St Paul Street, Suite 15, Rochester, NY 14605

P: 1.866.553.8536

V04302013