

## NEULOG PRESSURE SENSOR GUIDE



### NeuLog pressure logger sensor

#### NUL-210

The NeuLog pressure sensor can be used for any science experiment or activity which requires an accurate pressure measurement such as in the fields of Physics, Biology, Physiology, Microbiology, Botany, Weather Science, etc.

The pressure sensor comes pre-calibrated in the factory, and never requires recalibration.

Both the pressure and barometer sensors measure pressure however the pressure sensor is much easier to connect to lab ware such as syringes, beakers, and test tubes whereas the barometer sensor is ideal for testing atmospheric pressure in the open or in large containers. A hose with an attachment piece come connected to the sensor to make connections to various mediums easier.

Just a few of the thousands of possible experiments that can be done with the NUL-210 sensors are: Boyle's Law, Gay-Lussac's Law for ideal gases, weather phenomena, yeast and microbe fermentation, small animal exercise physiology, chemical reactions, and many more.

The pressure sensor can collect data using four different units of measure:

- Kilopascal (kPa): The SI unit of pressure.
- Atmospheres (atm): A pressure unit equal to one earth atmosphere at sea level (101.325 kPa)
- Pounds per square inch (psi): The unit of pressure which measures the amount of force (in pounds) a gas or liquid exerts across one square inch
- Bar: A non-SI unit equal to exactly 100 kPa

#### Quick start procedure:

##### PC or Mac Computer

##### Materials needed:

- NUL-210 Pressure Sensor
- USB-200 USB Module

- A USB to mini USB cable (which comes with the USB-200)

Your pressure sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via a USB to mini-USB cable. Please note that you cannot plug the pressure sensor directly into the computer.

Resident PC software and browser based software can be downloaded for free at [www.NeuLog.com/download](http://www.NeuLog.com/download) as well as a full software user guide. **Note:** Make sure not to download and install both types of software they will conflict on the computer.

##### Procedure:

1. Install the NeuLog software
2. Connect the USB-200 module to the PC or Mac
3. Connect the pressure sensor to the USB-200 module (they directly plug together). Please note there is no calibration required for this sensor.
4. Open the NeuLog software.
5. Once a pressure sensor logo appears on the left side of the screen the probe has been automatically identified and you can begin experimentation.
6. If the pressure sensor is not automatically identified then click the "Search for sensors" icon to find the sensor.
7. Select the "On-line experiment" button; this will open a graph below.
8. Click on the "Module setup" button located on the pressure sensor icon in the module window to change the sensor settings if need be.
9. Click on the experiment set up button to change the experiment settings if need be (experiment duration for example).
10. The pressure sensor will give a live reading in the box to the left of the screen while plugged in.
11. To run an experiment and collect data click "Run experiment".
12. To end data collection early, click "Stop experiment".

##### Tablet, smart phone device

##### Materials needed:

- NUL-210 Pressure Sensor
- WIFI-201 WIFI module
- BAT-200 Battery

## NEULOG PRESSURE SENSOR GUIDE



Your pressure sensor needs to be connected to a WIFI-201 module. The WIFI-201 module will create a closed NeuLog wifi network which will stream the NeuLog data to a device of your choosing. Once your device is wirelessly connected to the NeuLog network you can run experiments and collect data through a browser of your choosing.

### Procedure:

1. Connect the pressure sensor directly to the left side of a WIFI-201 module (no wires required).
2. Connect a BAT-200 module to the right side of the WIFI-201 module.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to micro USB charger (such as a typical cell phone charger). The WIFI-201 module will run for 15-75 minutes (depending on the sensor) without being plugged in.
4. For further WIFI-201 instructions or the WIFI-201 quick start guide please visit [www.NeuLog.com/download](http://www.NeuLog.com/download).
5. The WIFI-201 indicator lights will flash; take no action until the LED to the far left turns blue.
6. Take your tablet or smart phone and go to the Wi-Fi settings and select the NeuLog network which matches the WIFI module ID found on the back of the WIFI-201 device. (NeuLogXXXX for example)
7. Give your device 1-2 minutes to connect to the WIFI-201 network.
8. Once the device is connected go to your browser and type in the website [wifi201.com](http://wifi201.com) into the URL bar, then wait for 30-60 seconds.
9. You will see a "Control mode" icon in the browser, click on this icon.
10. The browser will then load a new screen and begin to auto detect the sensors, this can take a few minutes.
11. If the browser does not auto detect, select "Search for sensors"
12. Once the sensor is found you will see an icon on the left side of the screen for the pressure sensor. The icon will display data in real time.
13. To set the pressure sensor's settings click on "Module setup" button located on the pressure sensor icon on the left side of your screen.
14. To change the experimental settings click the "Experiment setup" button (experiment duration for example).
15. To run an experiment and collect data click "Run experiment".

### Operation with NeuLog viewer

#### Materials needed:

- NUL-210 Pressure Sensor
- VIEW-101 Viewing Module
- BAT-200 Battery

#### Procedure:

1. Connect the pressure sensor to the left side of the viewer.
2. Connect the BAT-200 module to the right side of the viewer.
3. The VIEW-101 will auto detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data click the run button (little green person).

### Off-line experiments

(Off-line experiments are for when you do not have a sensor connected directly to a computer, tablet, smartphone, or NeuLog viewer).

#### Materials needed:

- NUL-210 Pressure Sensor
- BAT-200 Battery

#### Materials needed to configure your offline experiment:

- USB-200 USB Module or WIFI-201 WIFI module or VIEW-101 Viewing Module
- A USB to mini USB cable (which comes with the USB-200)

#### Procedure:

1. Connect the pressure sensor directly to the left side of a charged BAT-200 module.

## NEULOG PRESSURE SENSOR GUIDE



2. When ready to collect data press the “Start/Stop” button on the pressure sensor, a red light will turn on for the duration of the experiment.
3. To change the experiment settings first connect to your chosen device (PC, Mac, tablet, smart device) and select the “Module setup” button. (For more information on how to change the experiment settings view quick start procedure section for your chosen device in this document.)
4. After the experiment has concluded (when the “Start/Stop” button is pressed again or the data collection period ends) connect the sensor to your chosen device. (full instructions in each device’s quick start procedure section)
5. Open NeuLog software.
6. Click the “Off-line experiment” button.
7. Click the “Load data from sensors” button.
8. Select which experimental data to upload. (5 experimental runs can be stored at one time on the NUL-210 pressure sensor)

### Restoring sensor’s factory default settings:

#### Procedure:

1. Connect the USB-200 to a PC, Mac, tablet, or smart device.
2. Connect the pressure sensor to a USB-200 module (they directly plug together). Please note there is no calibration required for this sensor.
3. Open the NeuLog software.
4. Click the “Tools” icon.
5. Click “Restore sensor’s factory defaults” icon.
6. If prompted to clear the graph, click either the “Off-line experiment” or “On-line experiment” (whichever you are using).
7. Click “Clear experiment results”.
8. Resume from step 4.

### Included with sensor:

- NeuLog NUL-210 Pressure Sensor Guide (this document).
- Attached pressure hose with end piece for easy connection.

### Sensor specifications

	atm	psi	kPa	bar
Range and operation modes	0 to 7	0 to 100	0 to 700	0 to 7
ADC resolution	13 bit			
Accuracy	±1%			
Resolution	0.01	0.1	0.1	0.01
Max sample rate (S/sec)	100			

### Sensor features:

- Fully digital data.
- Rugged plastic ergonomic case.
- Push button switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).
- Pre-calibrated internal sensing equipment.
- Hose with attachment piece connected directly to the sensor body for easy connection to various mediums.
- **Note:** NeuLog products are intended for educational use.

### Videos and experiment examples:

- Videos, literature and other probes can be found at [www.NeuLog.com](http://www.NeuLog.com)
- <http://www.NeuLog.com/Products/pressureloggerSensor.aspx>
- **Boyle’s Law** - <http://www.neulog.com/Portals/3/Users/Experiment C-30 Boyle's Law Ver 3.4.1.pdf>

### Technical background:

The philosophy behind NeuLog’s plug and play technology is based on each sensor’s ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (°C, °F, Lux, %, ppm, for example).

## NEULOG PRESSURE SENSOR GUIDE



The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

Inside the pressure sensor's plastic body there is a flexible membrane which bends with atmospheric pressure differences. On one side of the membrane there is a vacuum which is compared against the other side of the membrane, the atmosphere. An electrical charge is generated based on the amount of flexing in the membrane. Temperature is taken into account using internal circuitry during the pressure calculation as to normalize the data for any thermal fluctuations.

### Maintenance and storage:

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the pressure sensor's body.
- After use, gently wipe away any foreign material from the pressure sensor.
- Store in a box at room temperature out of direct sunlight.
- Do not blow directly into the sensor (No liquids should enter the sensor).

### Warranty:

We promise to deliver our sensor free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your sensor will be repaired or replaced.

Thank you for using NeuLog!



Flexible, simple, fast, forward thinking.

W: [www.neulog.com](http://www.neulog.com)

E: [info@neulog.com](mailto:info@neulog.com)

A: 850 St Paul Street, Suite 15, Rochester, NY 14605

P: 1.866.553.8536